

InstantAgarose™ TABLETS

Innovative Nucleic Acid Electrophoresis Workflow Essentials

HydraGene



InstantAgarose™ Protocol

Presoak tablet 2minutes in blow buffer volume before heating.

GEL%	1Tab	2Tab	3Tab
1%	50ml	100ml	150ml
1.2%	42ml	83ml	125ml
1.3%	38ml	77ml	115ml
1.5%	33ml	67ml	100ml
1.8%	28ml	56ml	83ml
2.0%	25ml	50ml	75ml

HyAgarose™ Green Agarose Product

HyAgarose™, Low EEO (LE) Agarose, Multi-purpose, only green agarose product in market, manufactured using advanced proprietary technology excludes the use of organic solvents. HyAgarose™ is a line of cleaner agarose products with significantly reduced environmental impact at improved cost efficiency while delivering same reliable high quality user experience.

ActGene HydraGene

Tel: (973) 978-6688 | Fax: (732)469-0333

E-mail: sales@actgene.com

Rethink Agarose

InstantAgarose™ TABLET

- Pre-weighted tablet, no weighing required
- Fast dissolving
- Made with high purity **HyAgarose™ Green LE Agarose**, multi-purpose
- Blister packaging allow for full portability and easy dispensing
- **<1 min. safe hands-on time**

Other Agarose POWDER

- Loose powder, need to be weighed
- Longer dissolving process
- Vary by manufacturer
- Bulky bottle packaging
- >20 min. hands-on time

Order Information

R9012TAB-200 InstantAgarose™ Tablet, 0.5gx200 tabs, blister convenient pack

Related product:

HydraGreen™ Safe DNA Dye, non-toxic, non-mutagenic formula of green safe alternative to Ethidium bromide

ACT-IDMG04 HydraGreen™ Safe DNA Dye, 20,000X in Water, 1ml

