



HydraGene, Co., Ltd.  
PO.Box 7011  
Piscataway, NJ 08855-7011  
Tel: 973-978-7888  
Web Site: [www.hydragene.com](http://www.hydragene.com)

China:  
E-mail [sales@sunmabio.com](mailto:sales@sunmabio.com)  
Tel: +86-592-6535930

## HydraGreen™ Safe DNA Dye, 10,000X in Water

### Product contents:

P/N: HDG-IDMG110

1x1ml

Sufficient for approximately 200x50ml or 100x100ml agarose gels

### Introduction:

HydraGreen™, 10,000X in water, is a non-carcinogenic and non-toxic alternative to Ethidium bromide used for the detection of nucleic acids in agarose gels. It is as sensitive as Ethidium bromide. There is also no toxic DMSO as HydraGreen™ is supplied in water.

HydraGreen™ has fluorescence excitation maxima at 295 nm and 490 nm. The fluorescence emission maxima is similar to EtBr when bound to DNA – at 530 nm.

### Protocol:

#### PRE-STAIN :

Add HydraGreen™ to melted agarose when the agarose has cooled to 50 to 60°C.

Use 8-12 µl per 100 ml of agarose or 4-6 µl per 50 ml of agarose

**IMPORTANT : HydraGreen™ is supplied in 10,000X concentration in water**

Or

#### POST-STAIN :

Use 20 to 30 µl per 100 ml of staining solution. Same as when using ethidium bromide. For an average gel thickness of about 7 mm, stain 30 minutes, followed by a destain of 30 minutes in water. Protect gel and staining solution from light with aluminum foil or place in dark.

#### DETECTION :

Detect bands under UV illumination (yellow or green gelatin- or cellophane filters is recommended for clearer bands) or non-UV LED illuminators such as Blue Light LED illumination.

### Shipping and Storage Conditions :

Shipped in RT

Store in RT for short term (1 month), Store in 4°C for long term. Protect from light.